

Compact Printer

AD-8121B



*Built for Electronic Balances, Counting Balances,
Platform Scales & Weighing Indicators.
Prints charts and offers a full range of Statistical
Functions on either AC or battery power.*



AND
A&D Company, Limited
www.aandd.co.jp

... Clearly a Better Value

Compact Printer

AD-8121B

Features

- A wealth of functions, including: Interval printing function, Statistical function, Calendar clock function and Chart print function
- Impact dot matrix printer for superior readability and long term durability
- A full range of statistical functions: Weight data, Total weight data, Counting data, Total counting data, Number of operations, Standard Deviation, Coefficient of Variant, Chart, Year, Month, Date, Hour, Minute and Second
- Changeable to dump mode while using statistical function without deletion of statistical function data
- Buffer memory of up to approximately 30 sets of data
- Interval printing with built-in Timer/Clock function: 5s, 10s, 30s, 1m, 5m, 10m, 30m
- Chart selectable to show change in weight, interval time, start time and stop time
- Compact size and light weight
- AC power supply and battery operation permits use anywhere
- 24-hour clock, calendar and pre-set functions(date and time) are stored in non-volatile memory (protected by a lithium battery)
- Accepts RS-232C and current loop input from A&D's Electronic Balances, Scales and Indicators
- Speedy printing

Specifications

Printer type ————— Dot impact type
 Character size ————— 5×7 dot, 2.5mm high, 16 characters/line
 Printing speed ————— Approximately 1 line/second
 Speed of paper feed ————— Approximately 1 line/second
 Print head life ————— About 500,000 lines
 Print paper ————— AX-PP143 (paper width 45mm, length 50m, diameter when rolled 65mm)
 Interface ————— RS-232C interface or current loop (2400bps)
 Power sources ————— AC adapter (provided as standard), or 4XAA alkaline batteries (battery life approximately 7000 lines)
 Dimensions ————— 180(W) × 160(D) × 80.5(H) mm
 Weight ————— About 400g (without print paper or batteries)
 Standard accessories — User manual, AC adapter, RS-232C connection cable (AX-KO-256A), print paper cover, printer paper axis, AA alkaline batteries (4)

Options

Current Loop Cable

Accessories (charged separately)

AX-PP143-S ————— Print paper (10 rolls)
 AX-ERC-05-S ————— Ink ribbon (5)
 AX-SW-128 ————— Foot switch
 AX-KO980-S100 ————— External input plug (necessary for foot switch connection)

[Note] When ordering this device, remember to specify whether RS-232C interface or a current loop cable will be used for connection and output.

Printing Examples

● Selectable printing modes

	Statistical Printing	Interval Printing	Chart Printing	
DATE 2004/03/17 TIME 10:00:21 WT 4032.58 g DATE 2004/03/17 TIME 10:00:37 WT 4036.27 g DATE 2004/03/17 TIME 10:01:46 WT 4029.15 g DATE 2004/03/17 QT 8251 PC DATE 2004/03/17 QT 7600 PC DATE 2004/03/17 QT 12389 PC TIME 13:10:25 WT 98.23 % TIME 13:11:43 WT 100.45 % TIME 13:14:37 WT 99.56 %	DATE 2004/03/17 TIME 15:00:00 ST. +00731.91 g ST. +00814.62 g ST. +05816.55 g ST. +00000.00 g ST. -00234.56 g WT 32.69 g WT 814.69 g WT 5816.55 g WT 0.00 g WT -439.68 g No. 34 WT +5836.94 g No. 35 WT +5816.55 g No. 36 WT +5746.28 g No. 37 WT +5721.61 g	No. 7 10:47:29 WT 9877.14 g No. 8 10:47:36 WT 9876.98 g No. 9 10:47:43 WT 9876.95 g No. 10 10:47:50 WT 9877.17 g No. 11 10:47:56 WT 9877.15 g DATE 2004/03/17 N 11 TOTAL 108647.66 g MAX 9877.48 g MIN 9876.64 g X 9977.06 g σ 0.235 g CV 0.01 % R 0.84 g	*INTERVAL TIME* 5 SEC. TIME 13:30:18 WT 9876.84 g TIME 13:30:23 WT 9876.92 g TIME 13:30:28 WT 9877.25 g TIME 13:30:33 WT 9877.00 g TIME 13:30:38 WT 9876.90 g TIME 13:30:43 WT 9876.59 g TIME 13:30:48 WT 9876.96 g TIME 13:30:53 WT 9877.19 g TIME 13:30:58 WT 9877.33 g *INTERVAL OFF*	*INTERVAL TIME* 5 SEC. DATE 2004/03/17 START 10:30:54 WT 2033.05 g 00h00m 00h04m STOP 10:35:47 WT 2033.57 g

Specifications are subject to change for improvement without notice.



...Clearly a Better Value

A&D Company, Limited

3-23-14 Higashi-Kebukuro, Toshima-ku, Tokyo 170-0013 JAPAN
 Telephone: [81] (3) 5391-6132 Fax: [81] (3) 5391-6148
www.aandd.co.jp

A&D ENGINEERING, INC.

1555, McCandless Drive, Milpitas, CA. 95035 U.S.A.
 Telephone: [1] (408) 263-5333 Fax: [1] (408) 263-0119

A&D MERCURY PTY. LTD.

32 Dew Street, Thebarton, South Australia 5031 AUSTRALIA
 Telephone: [61] (8) 8301-8100 Fax: [61] (8) 8352-7409

A&D INSTRUMENTS LTD.

Unit 24/26 Blacklands Way Abingdon Business Park,
 Abingdon, Oxon OX14 1DY United Kingdom
 Telephone: [44] (1235) 550420 Fax: [44] (1235) 550485

<German Sales Office>

Große Straße 13 b 22926 Ahrensburg, GERMANY
 Telephone: [49] (0) 4102 459230 Fax: [49] (0) 4102 459231

A&D KOREA Limited

Manhattan Bldg. 8F, 36-2 Yoido-dong, Youngdeungpo-gu, Seoul, KOREA
 Telephone: [82] (2) 780-4101 Fax: [82] (2) 782-4280